

iMatrix-411

Product No. AMS.892 041 350 µg
 Product No. AMS.892 042 1,050 µg

Version 001
 Store at 2-15 °C
 Protect from light

Background Information

Laminin-411 is well known to bind to the integrin α6β1 which is located on the cell surface. iMatrix-411 is recombinant Laminin411-E8 fragments which domain is known the binding region to integrin α6β1. iMatrix-411 increases the induction frequency to the blood vessel endothelial cells from pluripotent stem cells.

Content

Recombinant Human Laminin411-E8 Fragments

Amount

175 µg / tube (892 041: 2 tubes, 892 042: 6 tubes)

Form

Liquid solution

Product Information

iMatrix-411 is recombinant human Laminin411-E8 fragments expressed by CHO-S cell (Life Technologies).

Storage and Stability

The liquid solution is stable at +2 to +15 °C until the expiration date printed on the label.
 Protect from light.

iMatrix-411 is stable at 4 °C
 for 2 years from the manufacturing date.

Activity

The dissociation constant of the binding activity with integrin α6β1 is less than 10 nM.

Application

iMatrix-411 is able to be used as cell culture substrate for various cell types including ES/iPS cells.

Procedure

- 1) Dilute the solution with sterile PBS(-). Coat dishes with 0.5 µg/cm².
 * For example, for one well of a 6-well plate (9.6 cm²/well), add 9.6 µL of iMatrix-411 (4.8 µg) in 2 mL of PBS(-). Add 2 mL of diluted iMatrix-411 solution to the well.
- 2) Incubate for 1 h at 37 °C, 3 h at room temperature, or over night at 4 °C.
- 3) Remove remaining fluid from the coated surface. No rinse is needed.
- 4) Immediately plate the cells at desired density.
 * Don't allow the plate to dry.
 * The optimum coating concentration depends on cell lines, from 0.1 to 1.5 µg/cm².

References

Laminin 411 and 511 promote the cholangiocyte differentiation of human induced pluripotent stem cells
 Kazuo Takayama, et al.

Laminin-guided highly efficient endothelial commitment from human pluripotent stem cells
 Ryo Ohta, et al.

Regulatory Disclaimer

For life science research only. Not for use in diagnostic procedures.

AMSBIO | www.amsbio.com | info@amsbio.com

 **UK & Rest of the World**
 184 Park Drive, Milton Park
 Abingdon OX14 4SE, UK
 T: +44 (0)1235 828 200
 F: +44 (0) 1235 820 482

 **North America**
 1035 Cambridge Street,
 Cambridge, MA 02141
 T: +1 (617) 945-5033 or
 T: +1 (800) 987-0985
 F: +1 (617) 945-8218

 **Germany**
 Bockenheimer Landstr. 17/19
 60325 Frankfurt/Main
 T: +49 (0) 69 779099
 F: +49 (0) 69 13376880

 **Switzerland**
 Centro Nord-Sud 2E
 CH-6934 Bioggio-Lugano
 T: +41(0) 91 604 55 22
 F: +41(0) 91 605 17 85